


# A basic introduction to Dynamic Light Scattering (DLS) for particle size analysis

February 1st – 10:30 – 11:30 EST

## HOW TO LISTEN TO THE AUDIO

To listen to the webinar use your computer headphones or speakers

If you prefer to use your telephone to listen to the audio, please click on request audio in the participant panel

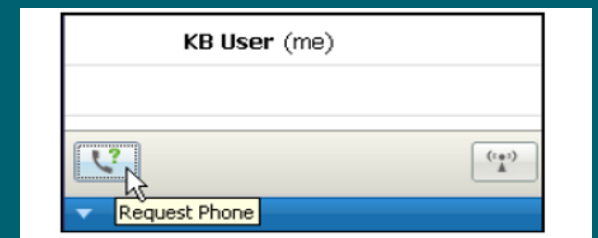



**Mike Kaszuba**

Technical Support Manager




**Craig Sagar**

Marketing Operations Manager



## DOWNLOAD A COPY OF TODAY'S PRESENTATION

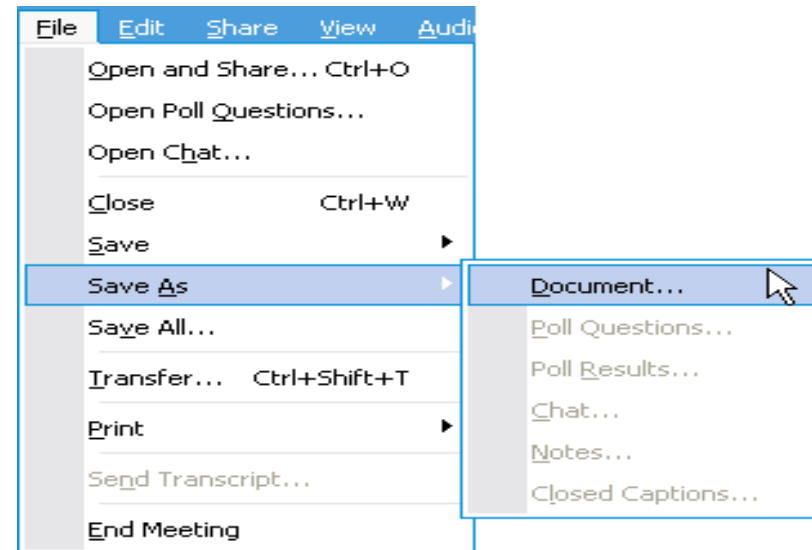

To save a copy of today's presentation, follow the instructions on the screen. Make sure you select \*.pdf in the "save as type" box.



**Malvern  
Panalytical**


# DYNAMIC LIGHT SCATTERING IN 30 MINUTES

## Particle and Molecular Size

Nano



$\mu$ V

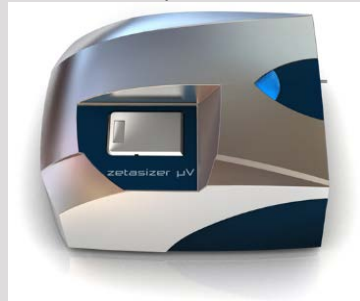

APS



**Dr Mike Kaszuba**

Technical Support Manager

E-mail: [michael.kaszuba@malvern.com](mailto:michael.kaszuba@malvern.com)

# Agenda



1. Brownian Motion
2. Correlation
3. Analysing the Correlation Function

# Agenda



1. **Brownian Motion**
2. Correlation
3. Analysing the Correlation Function

# Dynamic Light Scattering and Brownian Motion



- Non-invasive technique for measuring the size of particles and molecules in suspension
- Brownian motion is the **random** movement of particles due to the bombardment by the solvent molecules that surround them
- DLS measures the speed of particles undergoing Brownian motion
  - *Small particles diffuse rapidly*
  - *Large particles diffuse slowly*

# Brownian Motion



- Velocity of the Brownian motion is defined by the translational diffusion coefficient (D)
- The translational diffusion coefficient can be converted into a particle size using the **Stokes-Einstein** equation

$$d_H = \frac{kT}{3 \pi \eta D}$$

Where  $d_H$  = hydrodynamic diameter,  $k$  = Boltzmann's constant,  $T$  = absolute temperature,  $\eta$  = viscosity and  $D$  = diffusion coefficient



# Hydrodynamic Size




- Definition of Hydrodynamic Diameter ( $D_H$ ):


**The diameter of a hard sphere that diffuses at the same speed as the particle or molecule being measured**

- Dependent on
  - Ionic strength
  - Surface structure
  - Shape



# DLS Instrument Components

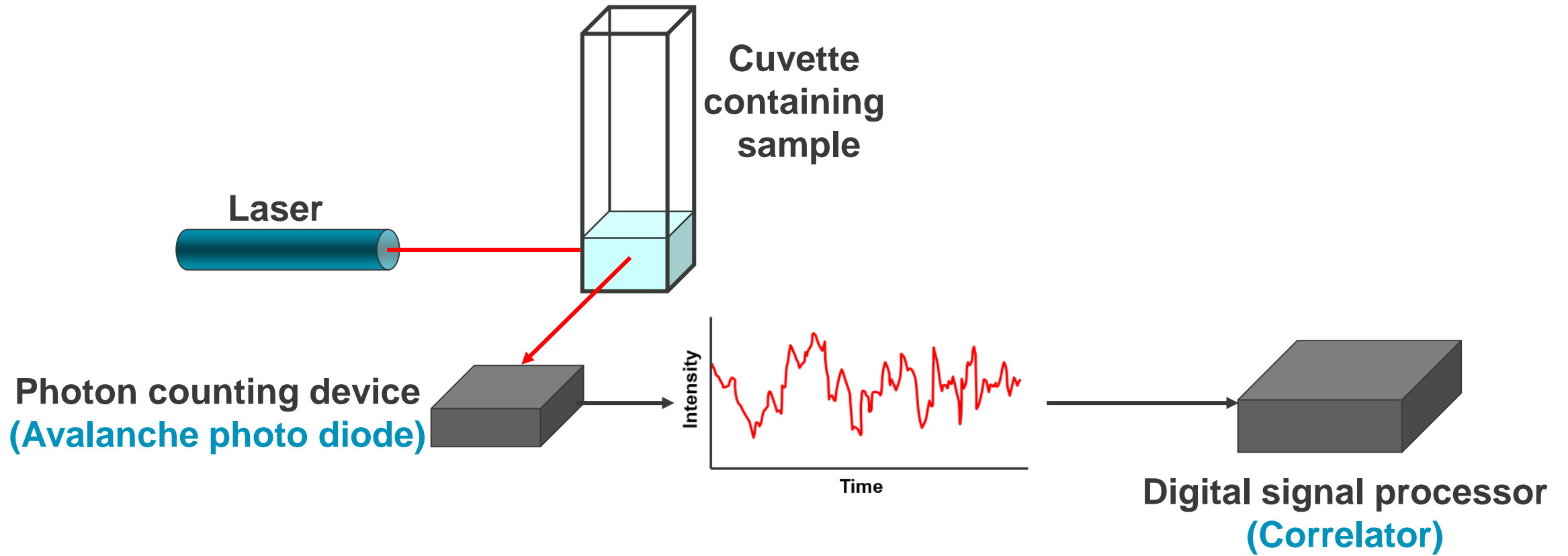

# Intensity Fluctuations and Brownian Motion

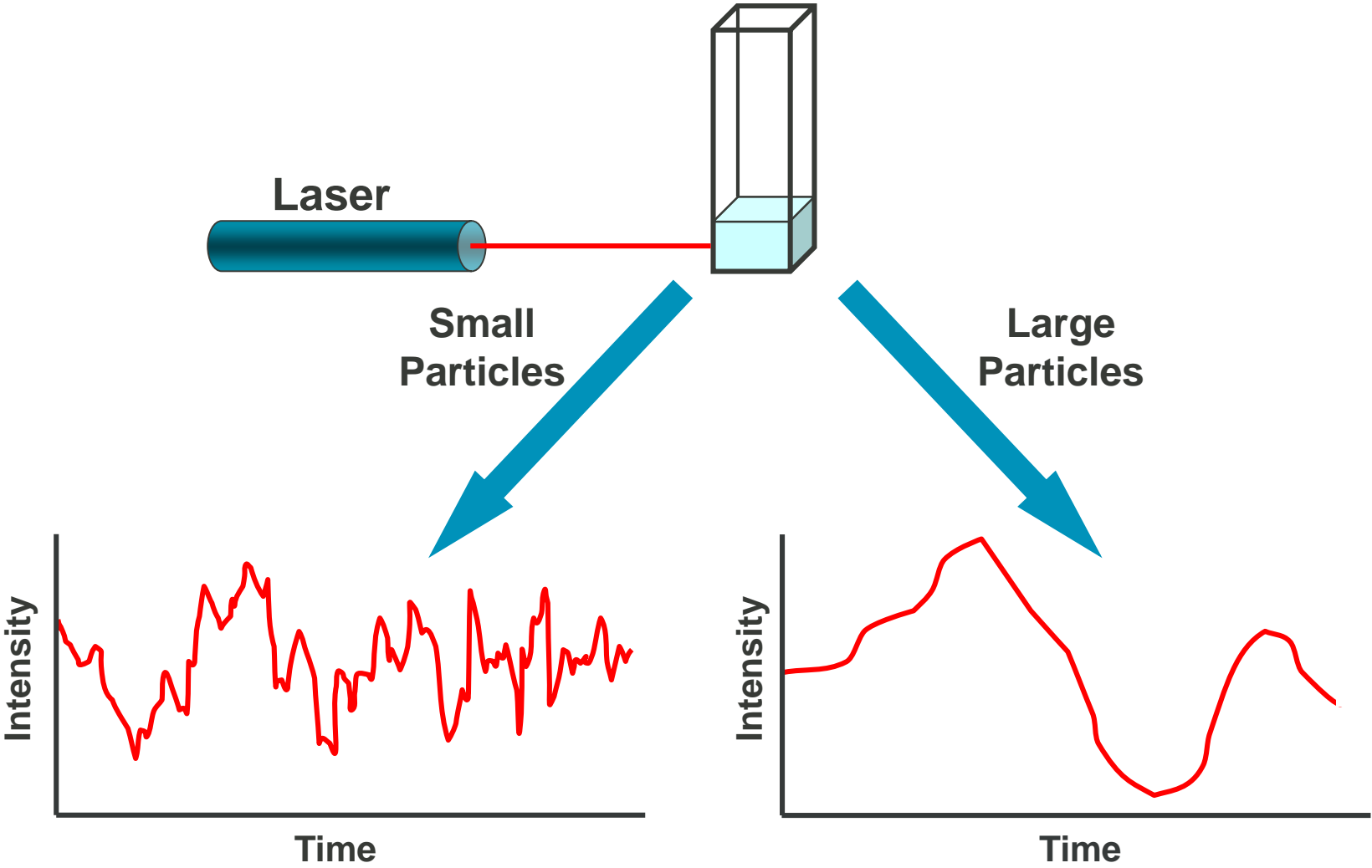

# Constructive and Destructive Interference

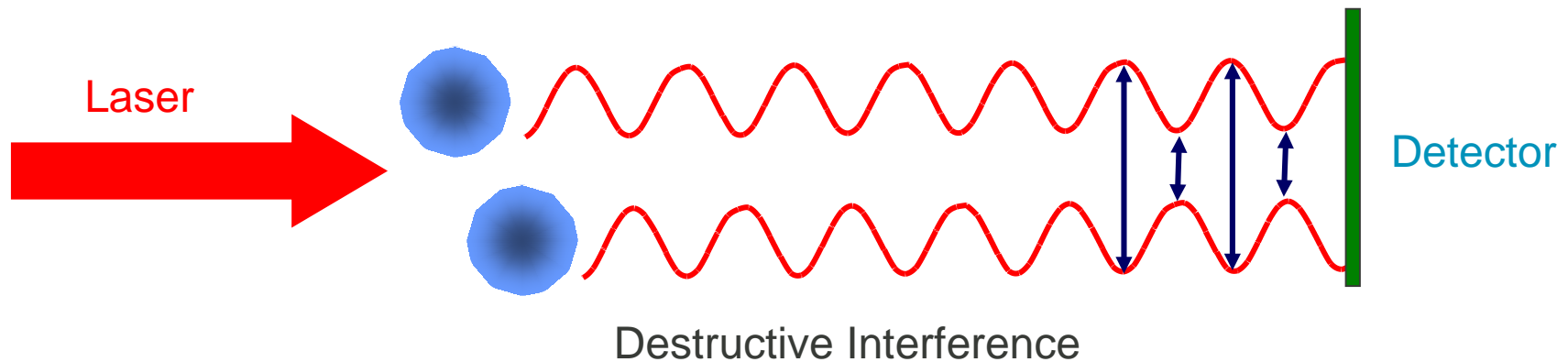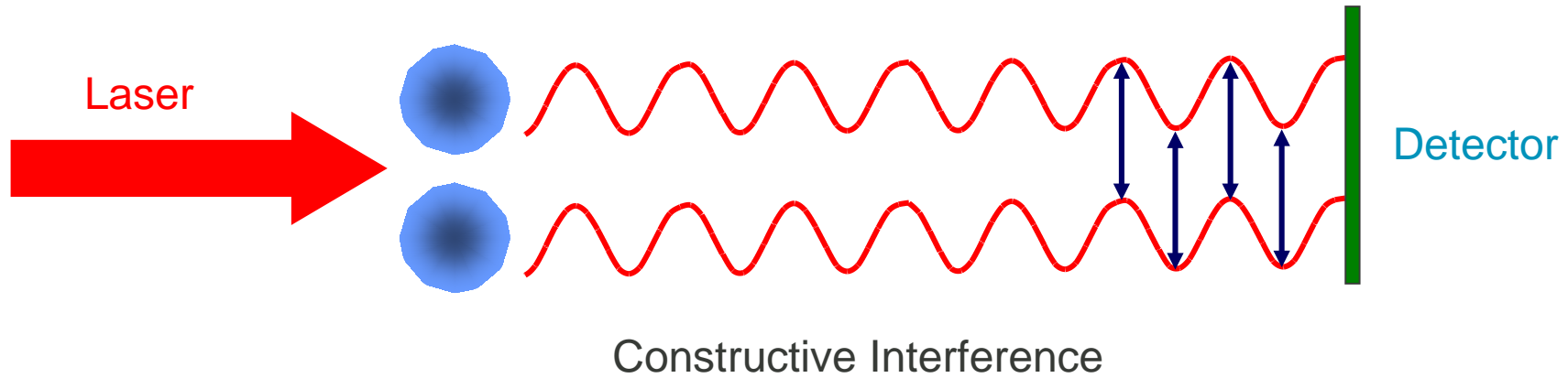

# Agenda



1. Brownian Motion
2. **Correlation**
3. Analysing the Correlation Function

# Correlation in Dynamic Light Scattering





- Technique for extracting the time dependence of a signal in the presence of “noise”
- Time analysis carried out with a correlator
- Constructs the time autocorrelation function  $G(\tau)$  of the scattered intensity according to





$$G(\tau) = \left\langle \frac{I(t_0) * I(t_0 + \tau)}{I(t_\infty)^2} \right\rangle$$

where  $I$  = intensity,  $t$  is the time and  $\tau$  = the delay time

# Correlation



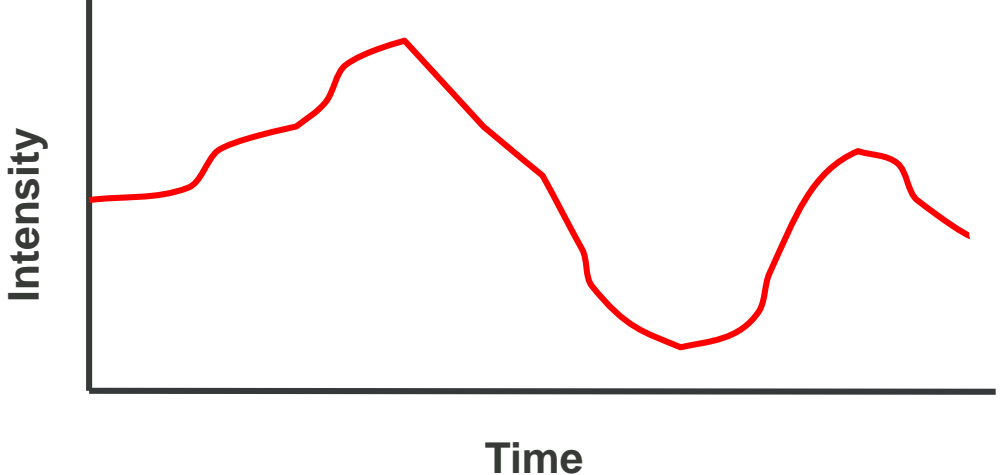
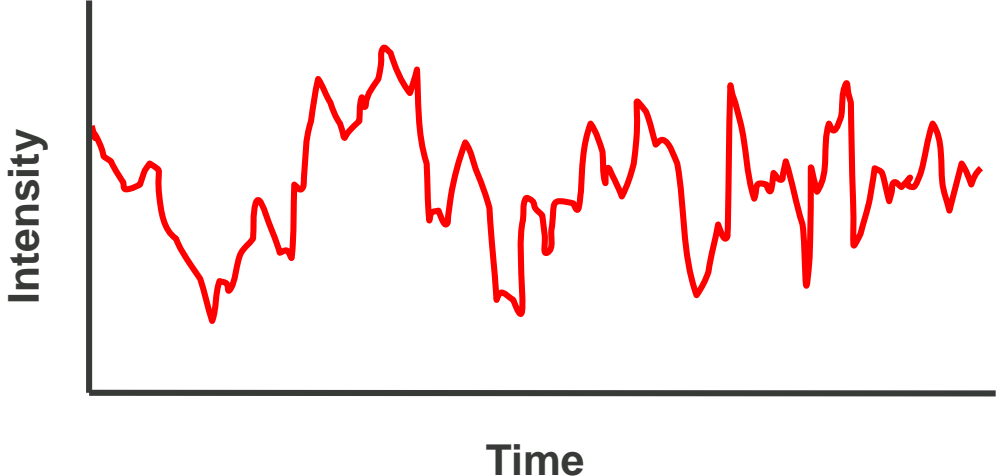


# Correlation

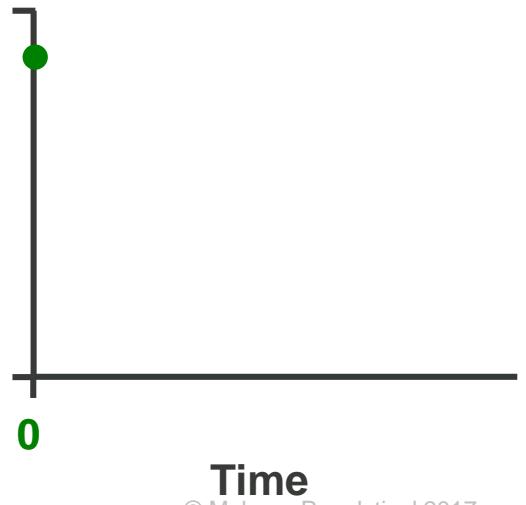
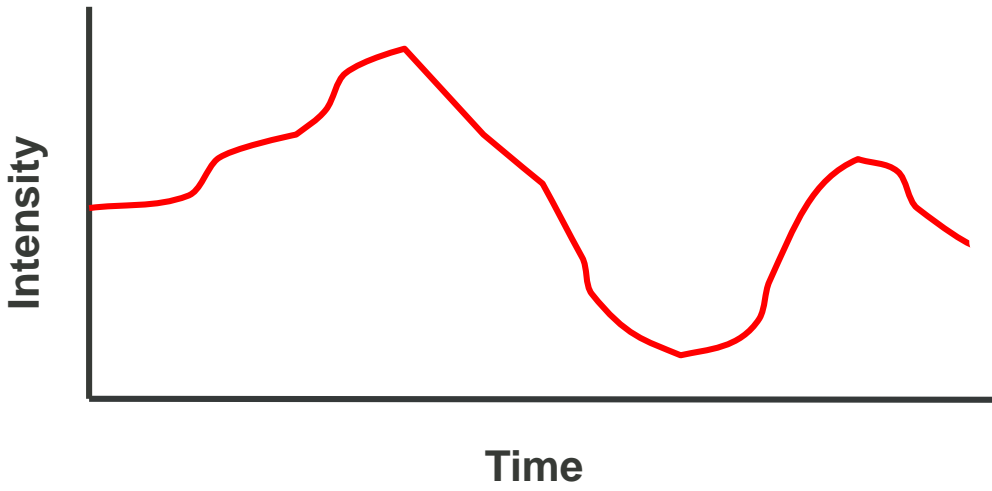
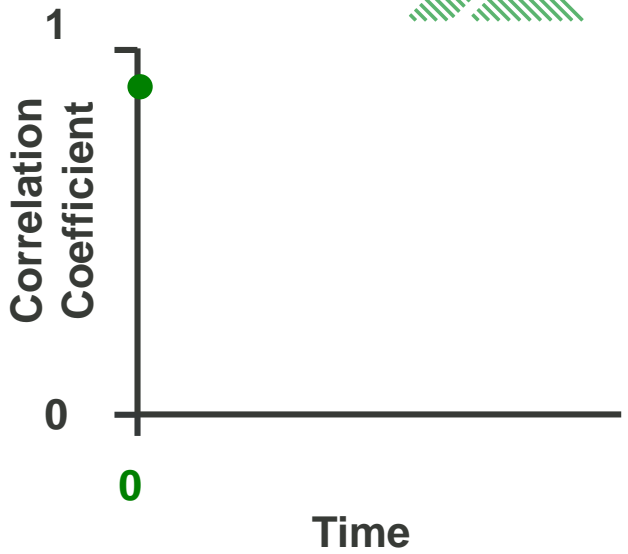
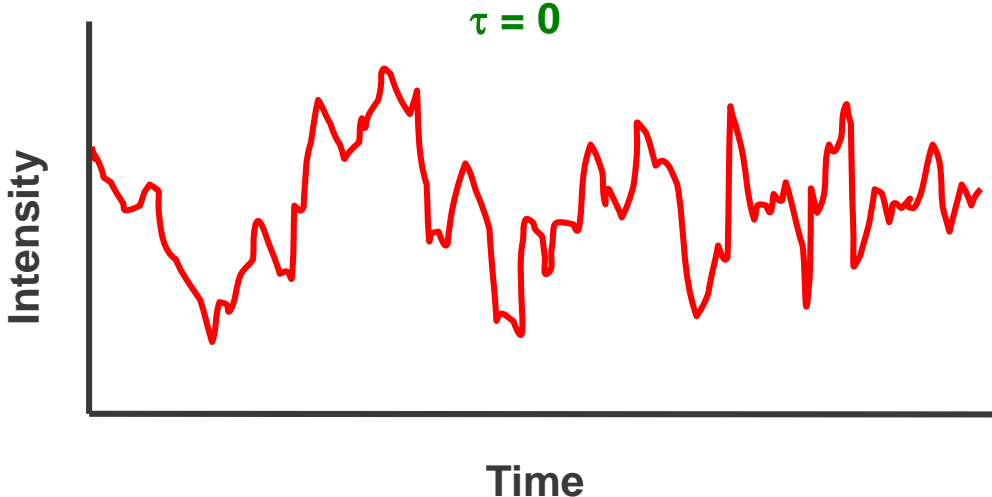








# Correlation

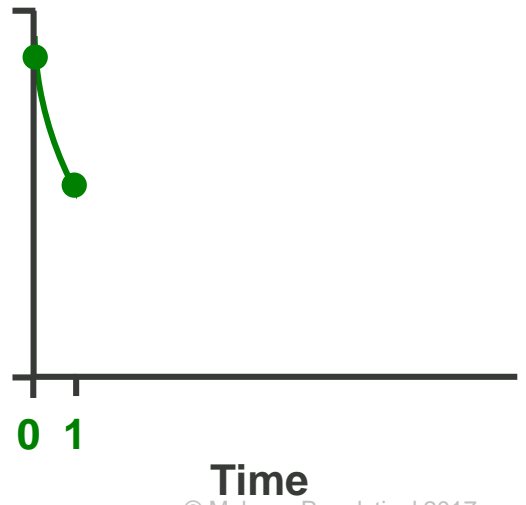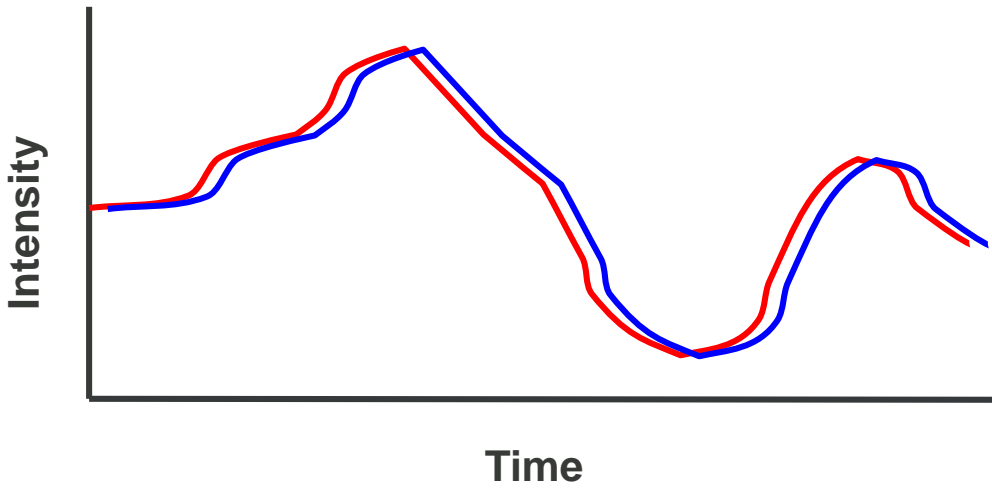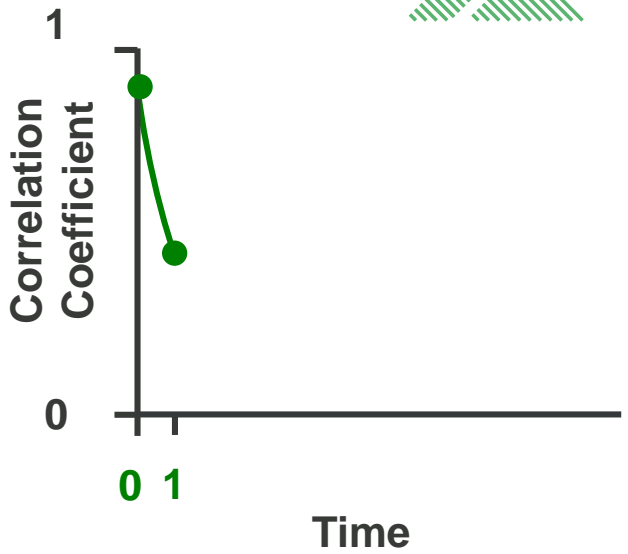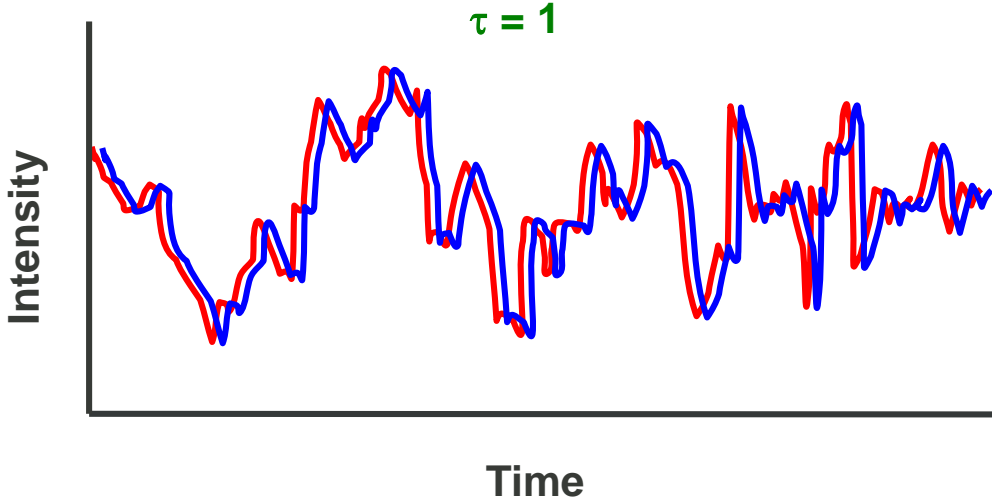





# Correlation

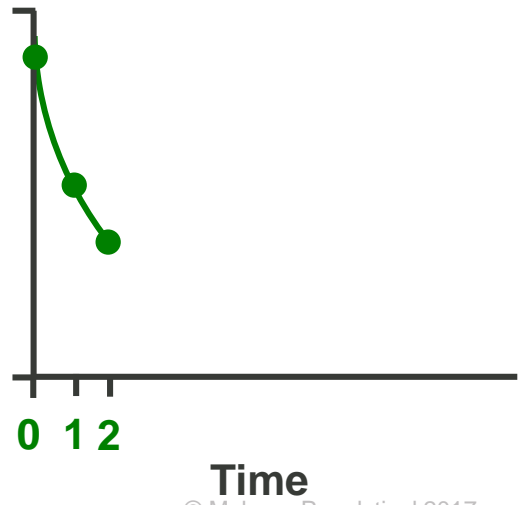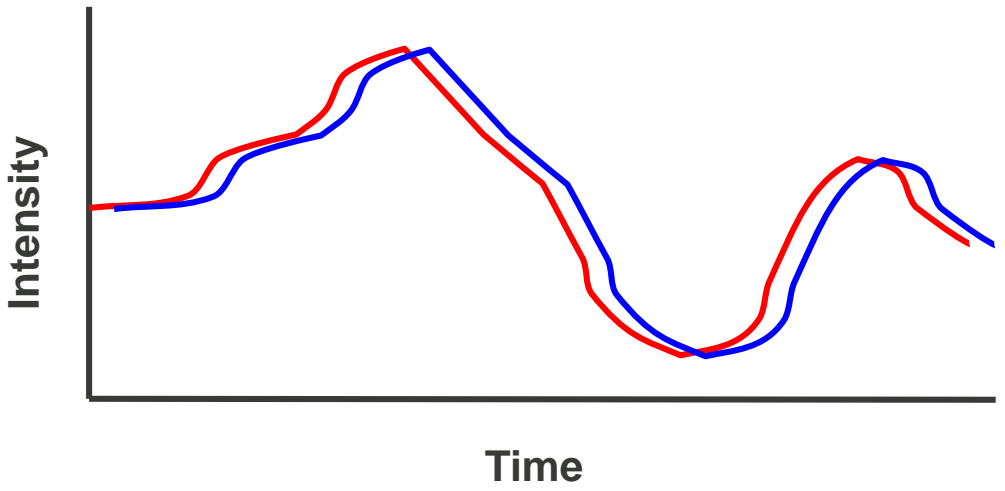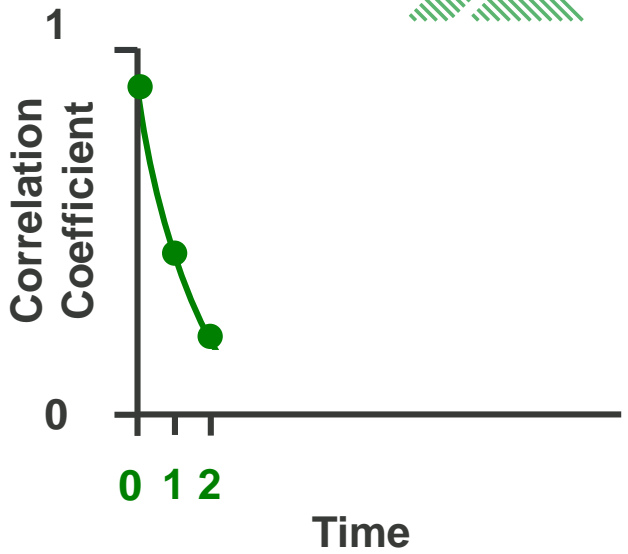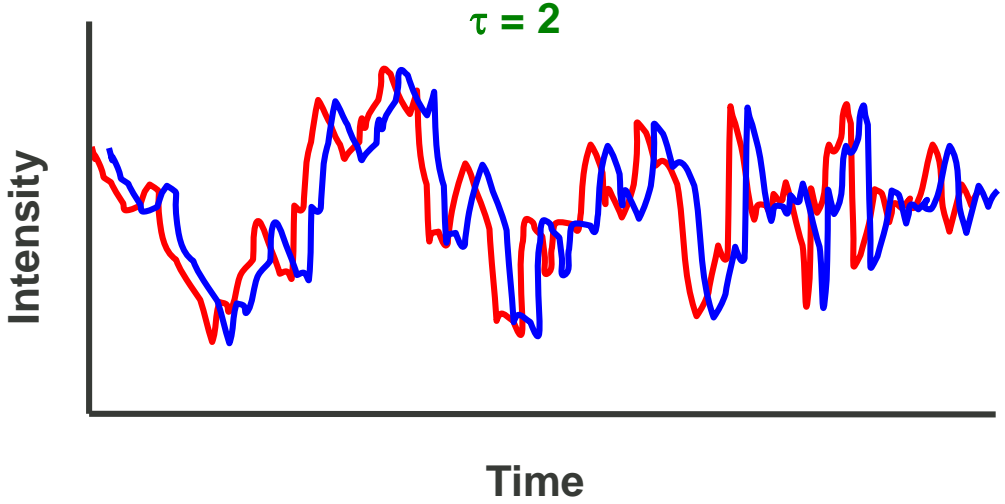





# Correlation

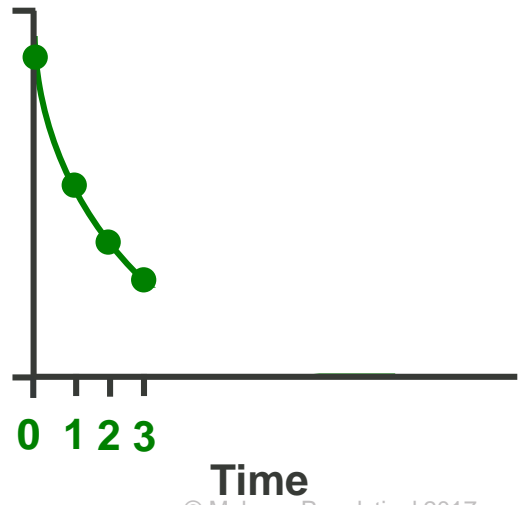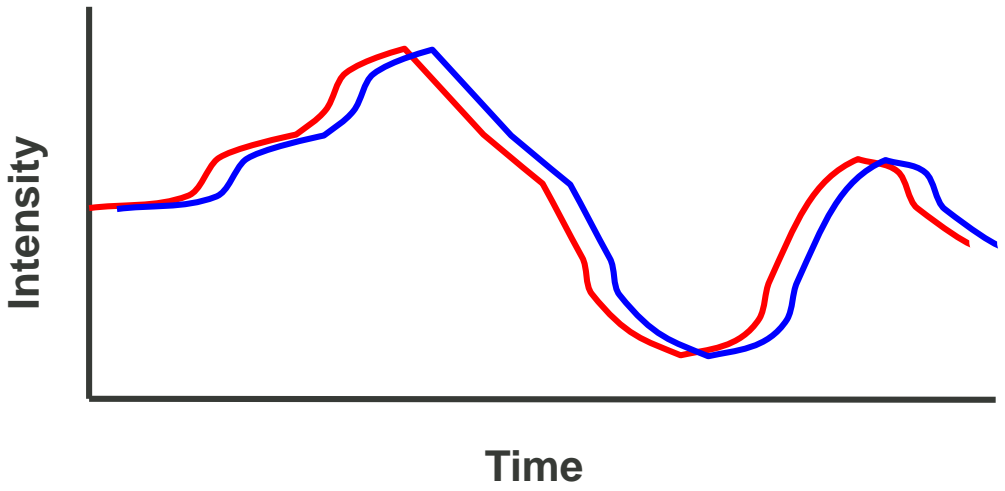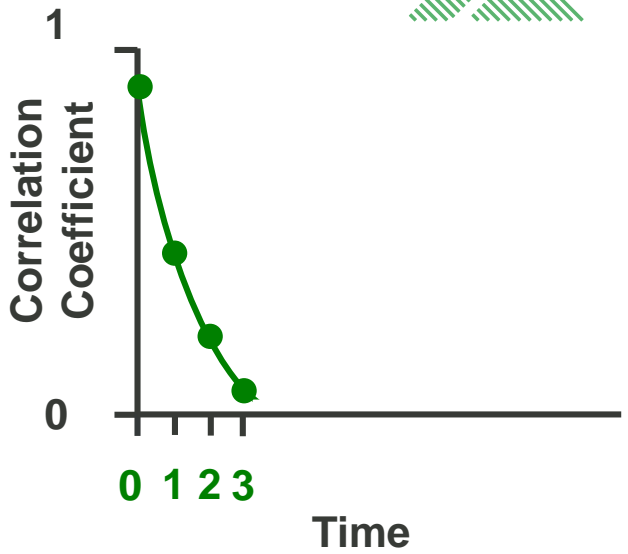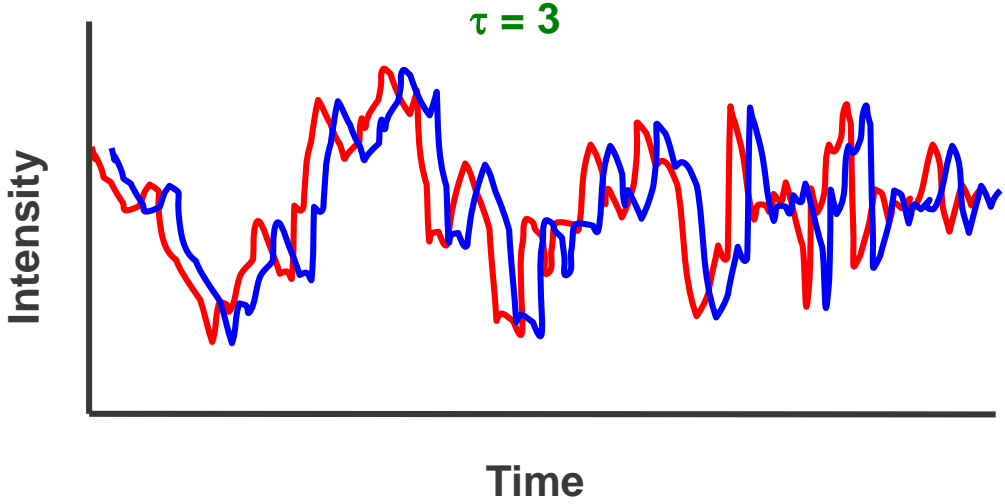


# Correlation

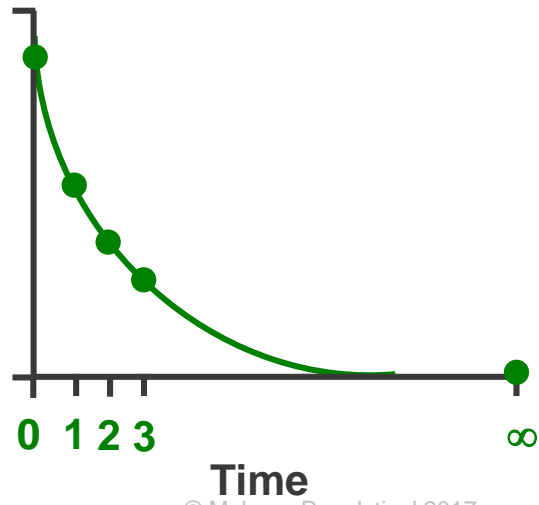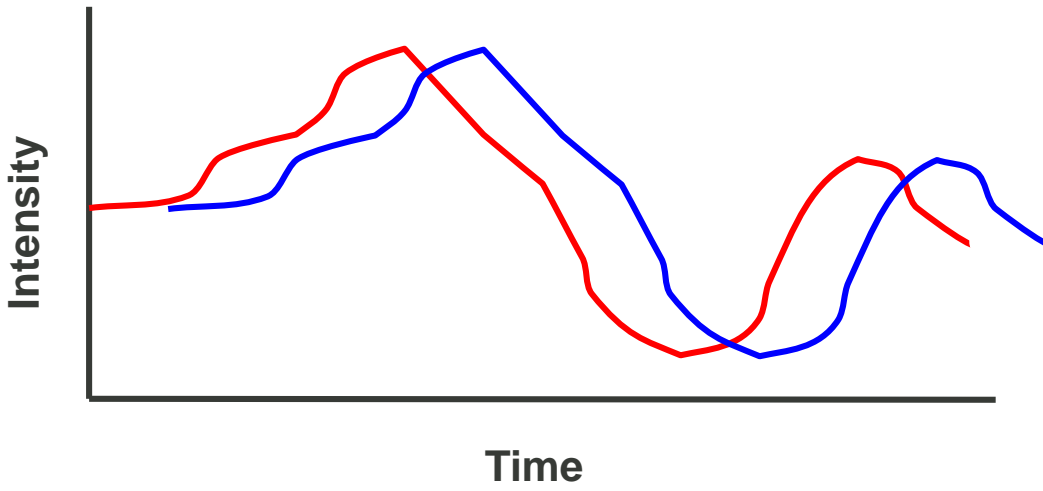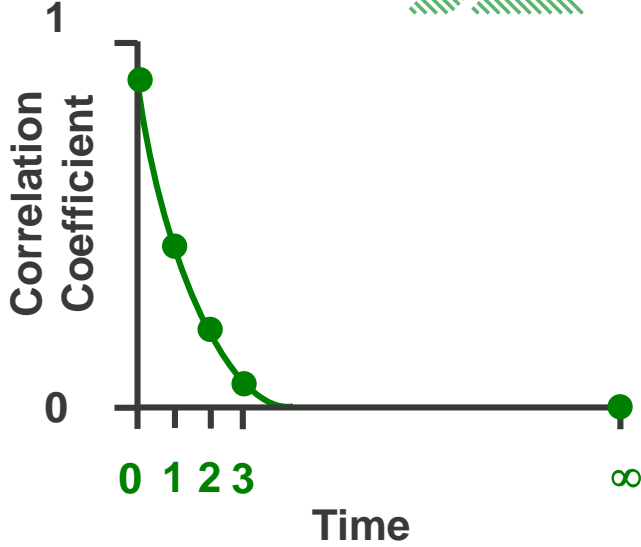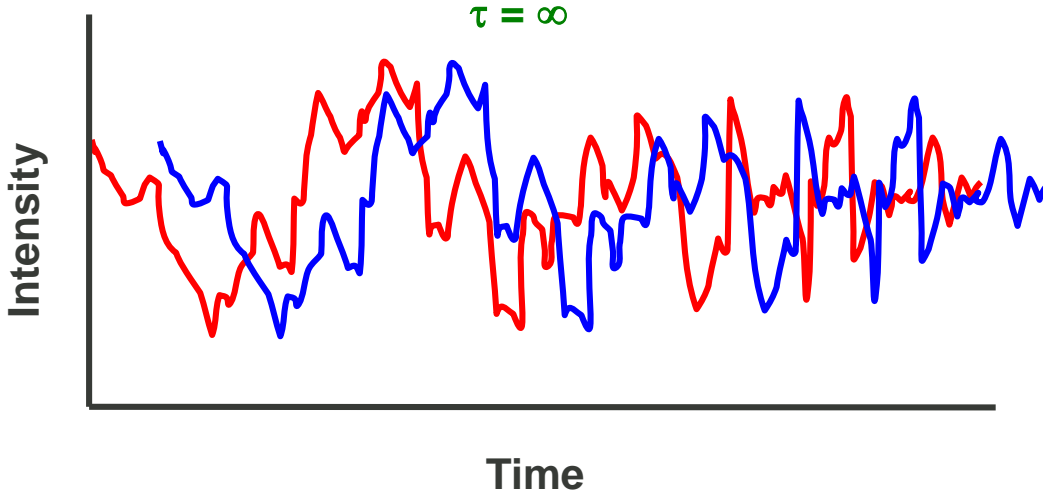

# Correlation

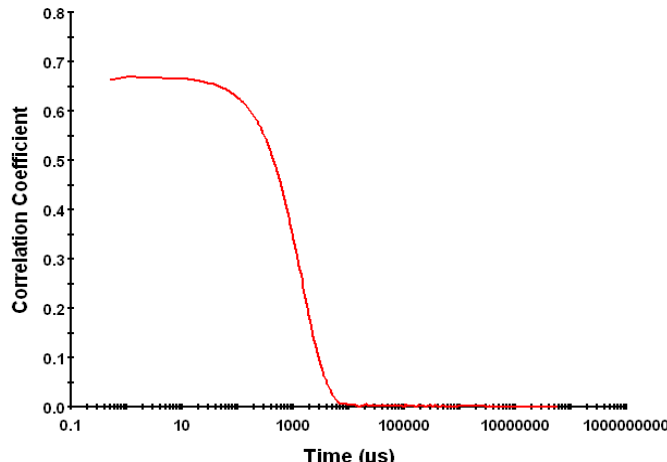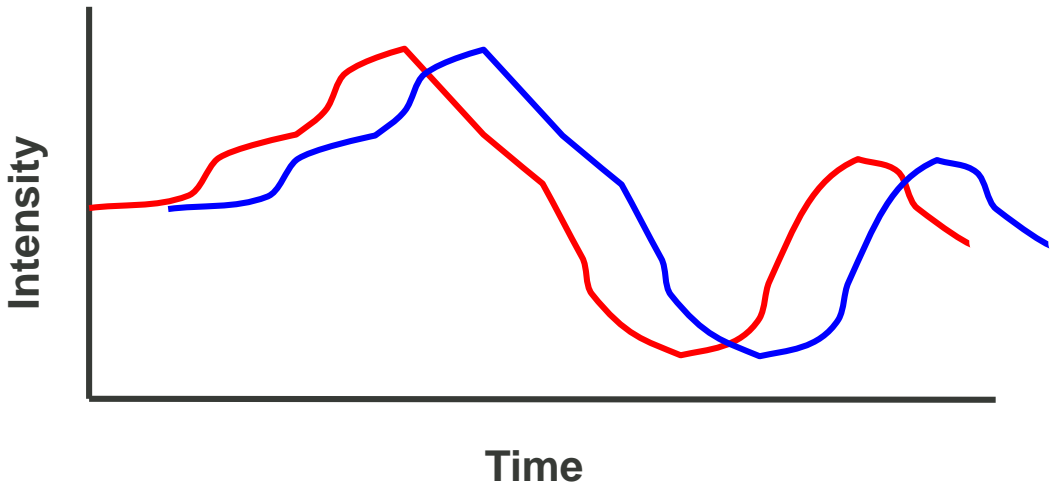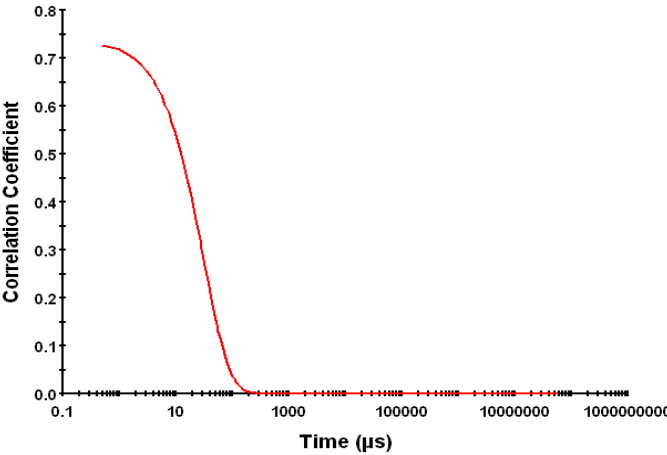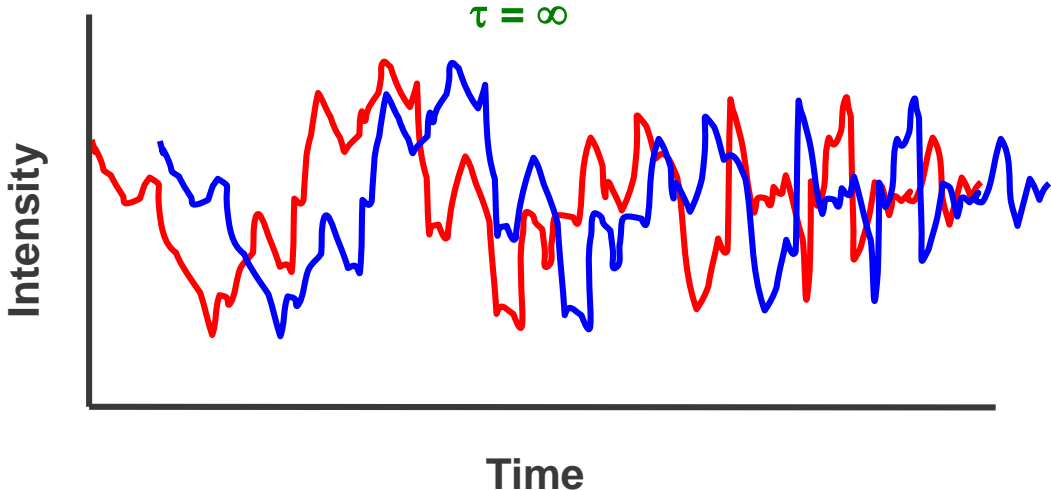

# Correlation Functions

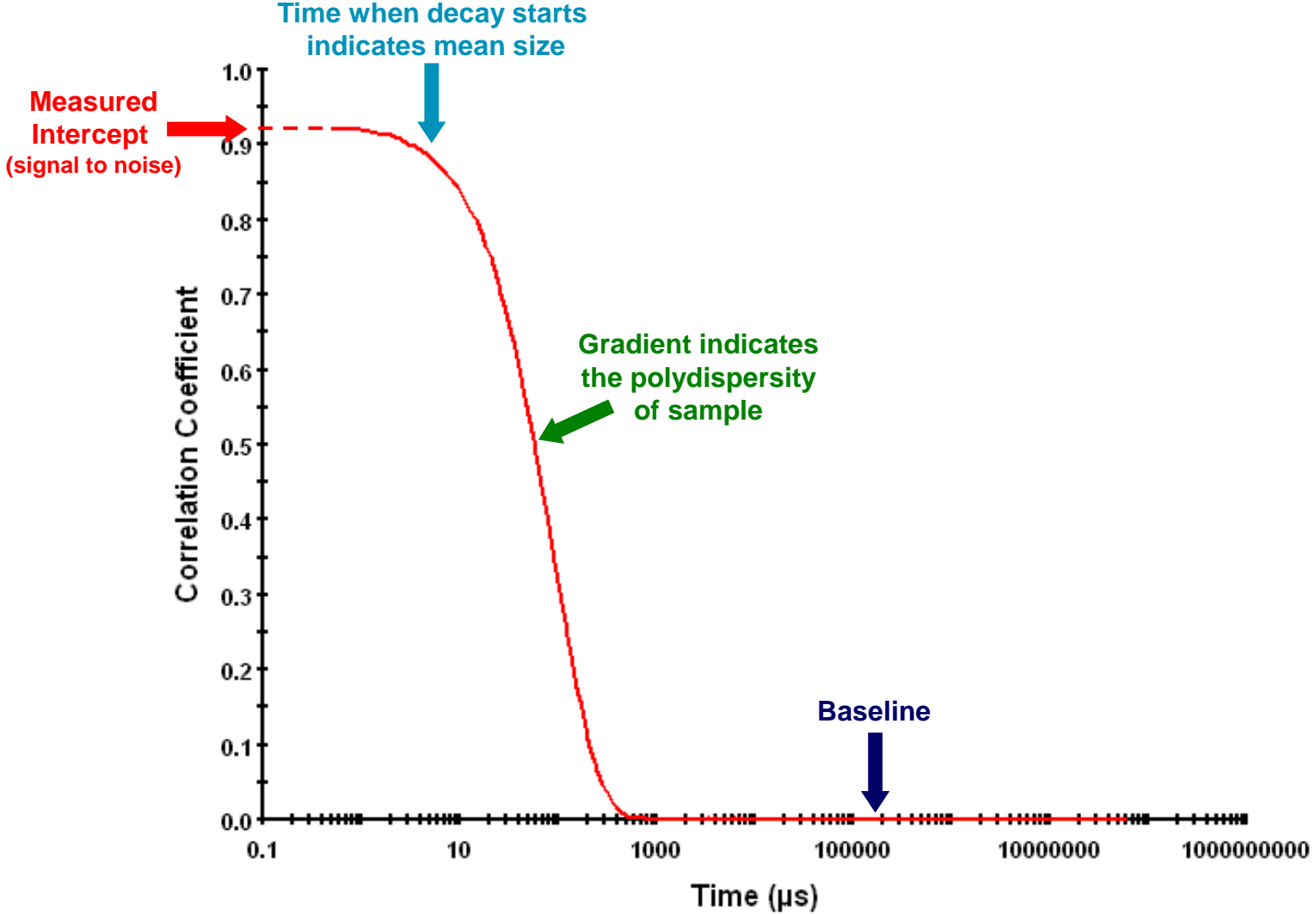

# Agenda



1. Brownian Motion
2. Correlation
3. **Analysing the Correlation Function**

# Correlation Functions



- The correlation function can be modelled with an exponential expression such as:

$$G(\tau) = B + A \sum e^{-2q^2 D \tau}$$

Where

B = baseline at infinite time

A = amplitude (or intercept)

q = scattering vector =  $(4\pi n/\lambda_0) \sin(\theta/2)$

n = dispersant refractive index

$\lambda_0$  = laser wavelength

$\theta$  = detection angle

D = diffusion coefficient

$\tau$  = correlator delay time



# Analysing The Correlation Function



- Correlation function contains the **diffusion coefficient** information required to be entered into the Stokes-Einstein equation
- The **diffusion coefficients** are obtained by fitting the correlation function with a suitable **algorithm**

# Analysing The Correlation Function




Two different analyses are performed:

## Cumulants analysis

- Mean size (z-average)
- Polydispersity index

## Distribution analysis

- Distribution of sizes

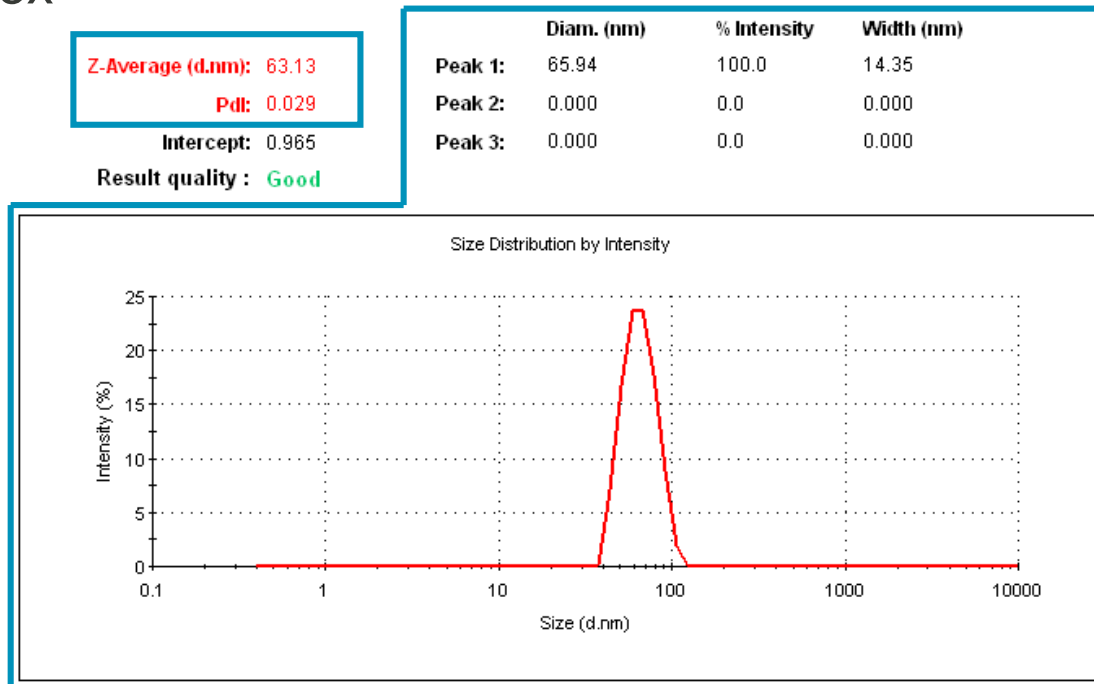

# Cumulants Analysis



- Defined in the International Standards ISO13321 (1996), ISO22412 (2008) and **ISO22412 (2017)**
- Only gives a **mean particle size (z-average)** and an **estimate of the width of the distribution (polydispersity index)**
- Only the dispersant refractive index and viscosity are required for this analysis

# The z-Average Diameter



- Definition of the z-Average Diameter ( $Z_D$ ):

**The intensity-weighted mean diameter derived from the cumulants analysis**

- Specific to light scattering
- Very sensitive to the presence of aggregates or large contaminants due to the inherent intensity weighting

# Polydispersity Index



- Definition of the Polydispersity Index (Pdl):

**A dimensionless measure of the broadness of the size distribution calculated from the cumulants analysis**

- Ranges from 0 to 1 in the Zetasizer software
- Values  $> 1$  indicate that the distribution is so polydisperse, the sample may not be suitable for measurement by DLS

# Zetasizer Distribution Algorithms



- **General Purpose**

- Suitable for the majority of samples where no knowledge of the distribution is available
- Will give broad, smooth distributions

- **Multiple Narrow Modes**

- Suitable for samples suspected to contain discrete populations
- Will give narrow peaks


- **Protein Analysis**

- Best suited for protein samples – will give narrow peaks – can be used for any sample type
- Automatically picks the optimal distribution

# Size Distributions in the Zetasizer Software




- Primary size distribution is **intensity**-weighted
- A plot of the **relative intensity of light scattered** by particles (on the Y axis) versus various **size classes** (on the X axis) which are logarithmically spaced

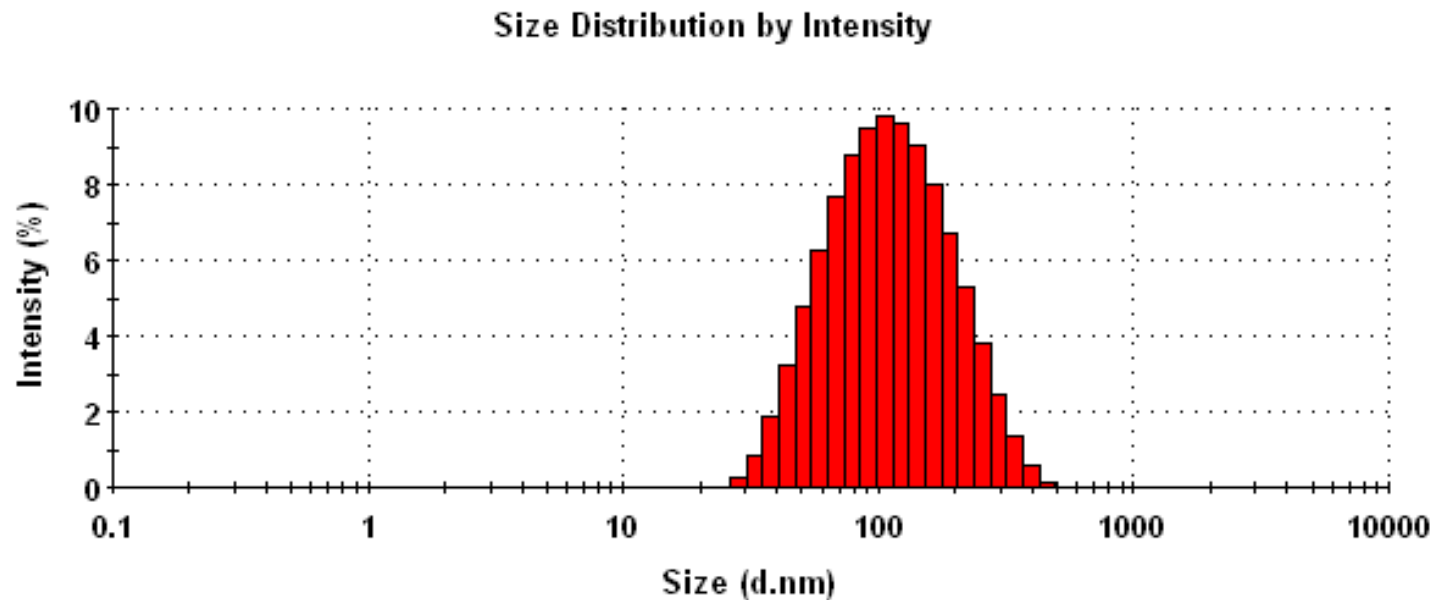

# Intensity Size Distributions



- Primary result
- Based upon the intensity of light scattered by particles
- Sensitive to the presence of large particles/aggregates /dust
- Only the dispersant viscosity and refractive index values are required

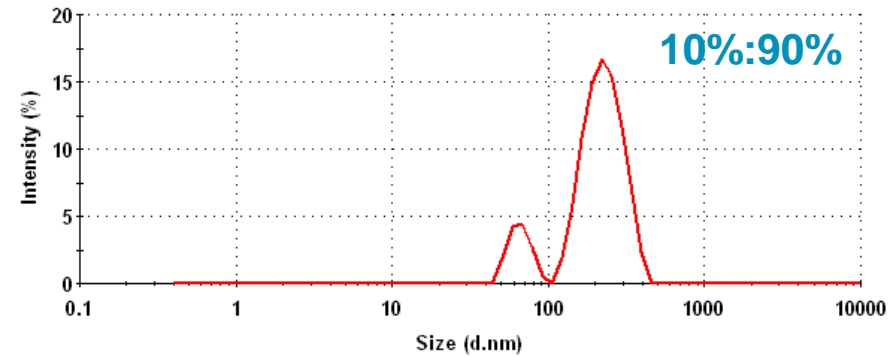






# Volume Size Distributions

- Derived from the intensity distribution using Mie theory
- Equivalent to the mass or weight distribution
- Particle optical properties required to make this transformation
  - Particle refractive index
  - Particle absorption

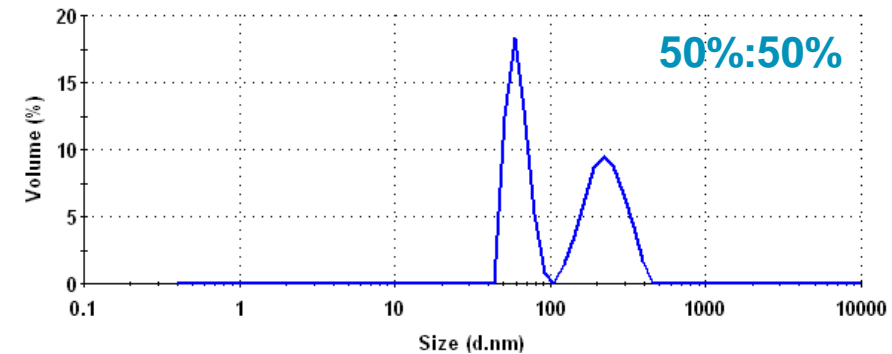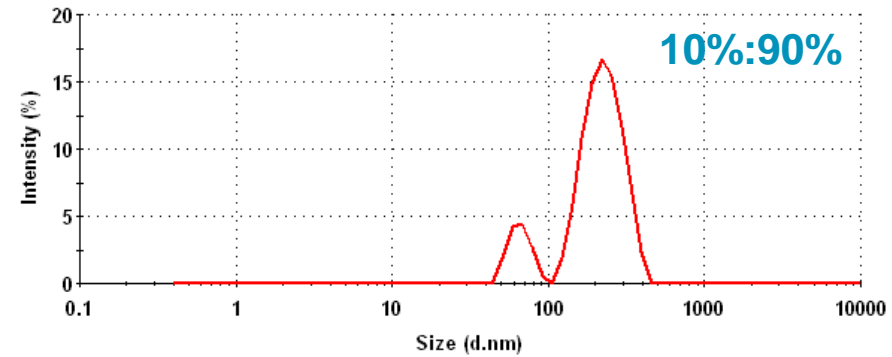

# Number Size Distributions

- Derived from the intensity distribution using Mie theory
- Particle optical properties required to make this transformation
  - Particle refractive index
  - Particle absorption

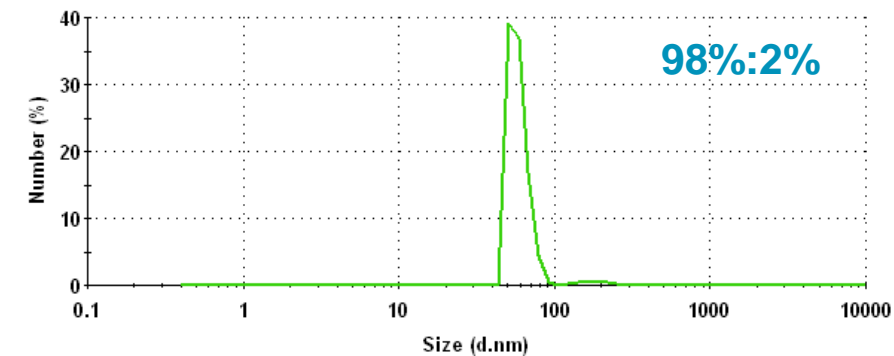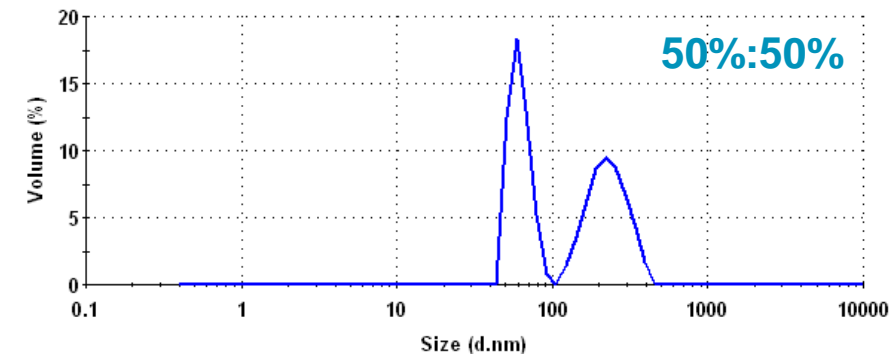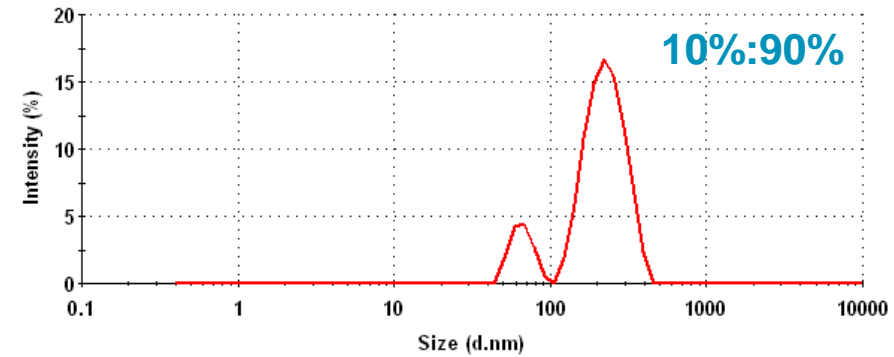

# Size Distributions From DLS



- DLS technique tends to overestimate the width of the peaks in the distribution
- This effect is magnified in the transformations to volume and number
- The volume and number size distributions should only be used for estimating the **relative amounts** of material in separate peaks as the means and particularly the widths are less reliable


# Volume/Number Distributions: Recommended Use



- Use the **Intensity PSD** for reporting the size of each peak in the distribution
- Use the **Volume** or **Number PSD** for reporting the relative amounts of each peak in the distribution

(Modal Size Report)

<b>Z-Average (d.nm):</b> 175.1	<b>Diam. (nm)</b>	<b>Width (nm)</b>	<b>% Intensity : Volume : Number</b>		
<b>PdI:</b> 0.191	<b>Peak 1:</b> 222.9	51.67	90.8	52.3	2.0
<b>Intercept:</b> 0.866	<b>Peak 2:</b> 57.07	7.544	9.2	47.7	98.0
<b>Result quality:</b> Good	<b>Peak 3:</b> 0.000	0.000	0.0	0.0	0.0

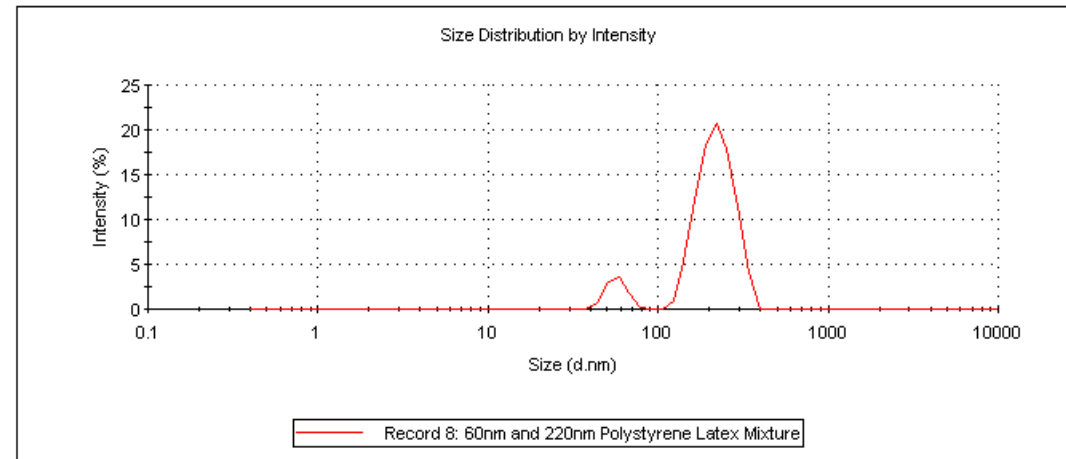



# Any Questions?

[michael.kaszuba@malvernpanalytical.com](mailto:michael.kaszuba@malvernpanalytical.com)

[craig.sagar@malvernpanalytical.com](mailto:craig.sagar@malvernpanalytical.com)